



# Risperidone-loaded PLGA sustained release implants

---

from benchtop to commercial scale

Patrick Duffy, Ph.D.

CRS 2024 Annual Meeting

July 8 – 12 Bologna, Italy

# collaborators and case study

---


# viatel™ polymers for implants - case study

#	Polymer type/family	Product Code	Batch	L : G (mol%)	Molecular Weight Mn, Mw (kDa)	IV (dl/g)	End-group
1	Poly D,L-lactide-co-glycolide (PLGA)	DLG 5002 E	2591607	49.4 : 50.6	9.0, 18.3	0.20	Ester
2	Poly D,L-lactide-co-glycolide (PLGA)	DLG 5002 A	2588968	49.4 : 50.6	8.8, 19.2	0.20	Acid
3	Poly D,L-lactide-co-glycolide (PLGA)	DLG 5005 A	2567373	49.7 : 50.3	29.8, 54.3	0.49	Acid
4	Poly D,L-lactide-co-glycolide (PLGA)	DLG 7502 A	2597443	73.9 : 26.1	7.8, 16.2	0.20	Acid
5	Poly D,L-lactide-co-glycolide (PLGA)	DLG 7509 E	2573112	74.5 : 25.5	61.9, 104.2	0.82	Ester
6	Poly D,L-lactide-co-glycolide (PLGA)	DLG 8507 E	2588392	84.5 : 15.5	49.9, 90.7	0.70	Ester
7	Poly D,L-lactide (PDLLA)	DL 03 A	2550133	100 : 0	28.5, 50.8	0.39	Acid
8	Poly D,L-lactide (PDLLA)	DL 09 E	2596557	100 : 0	57.6, 89.6	0.72	Ester



8 distinct grades of amorphous polyesters utilised for screening  
varying: L:G ratios, Mw / IV, end-group

# screening using VCM tool – no drug

checking the influence of melt processing on various viatel™ polymers

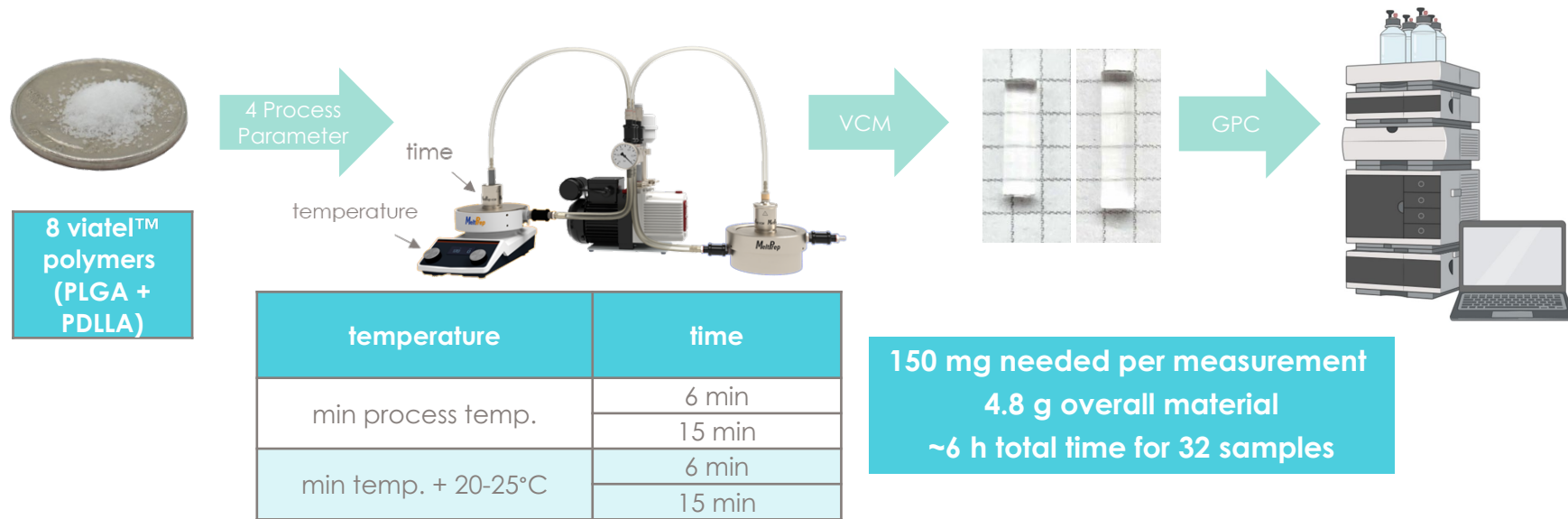


# process influence on viatel™ polymers

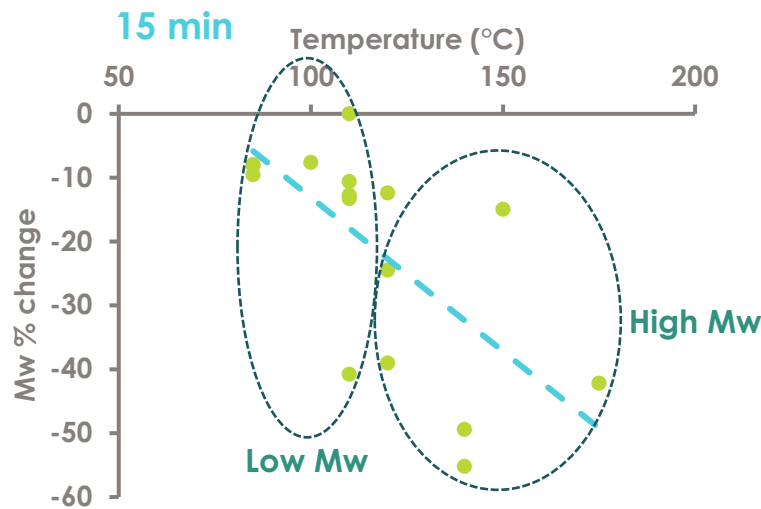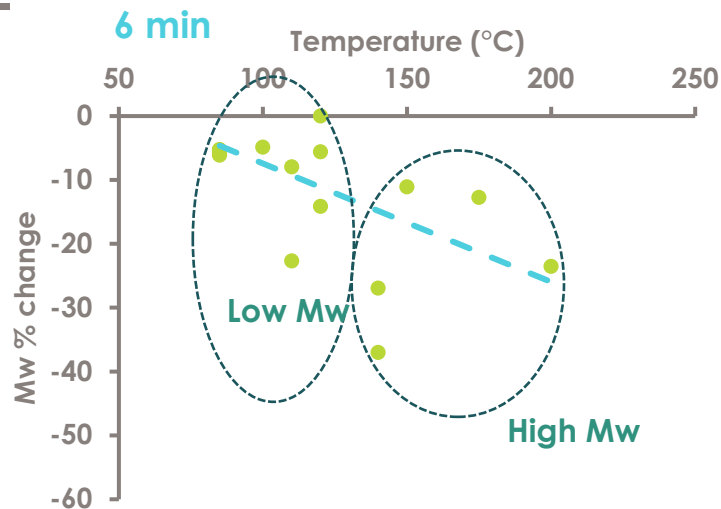

melt processing should be done at mildest conditions possible  
to minimize Mw decay



# process influence on viatel™ polymers

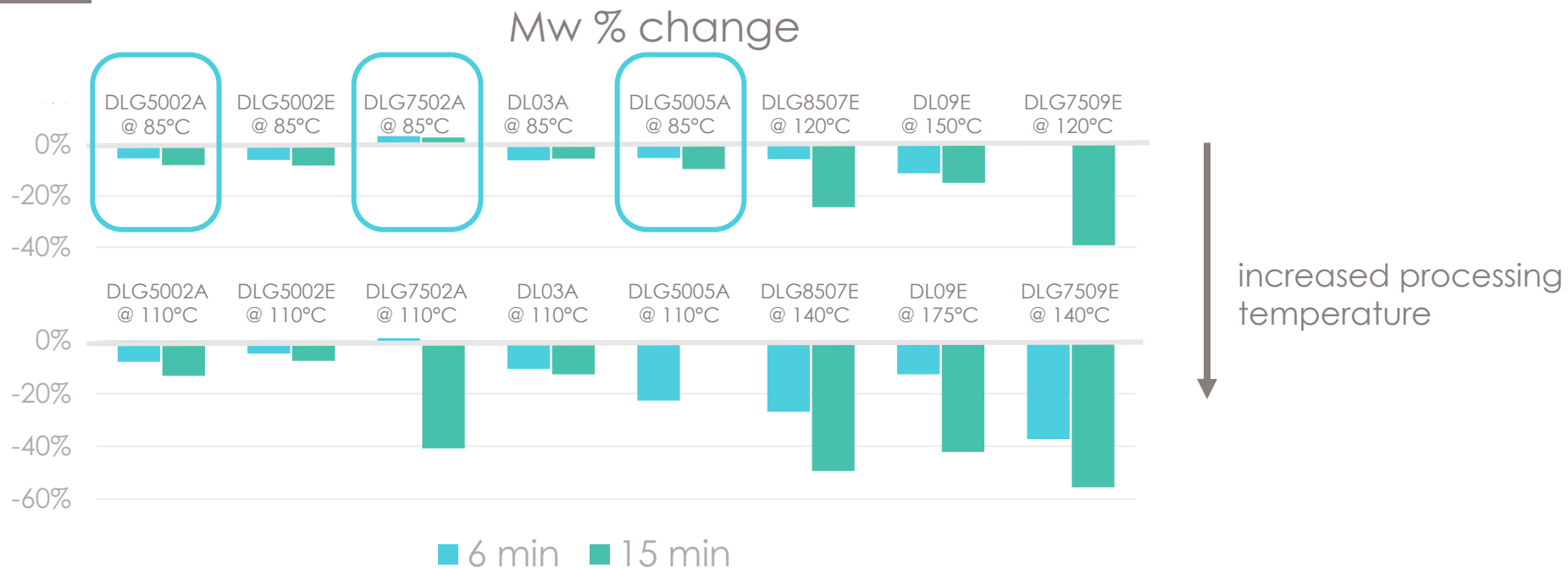

# implant fabrication using VCM tool

viatel™ PLGA

risperidone

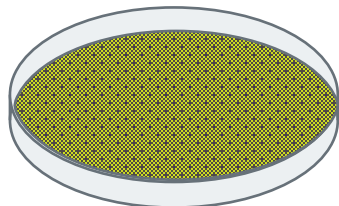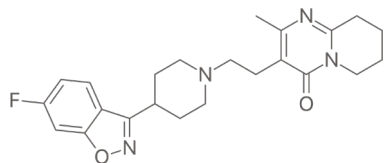

viatel™ PLGA + risperidone film



mild conditions  
100°C / 5 min




∅ 3 mm



viatel™ PLGA / risperidone implants

# implant characterisation – SEM & Raman

- SEM shows uniform, well-formed implants
- Raman shows good drug distribution throughout

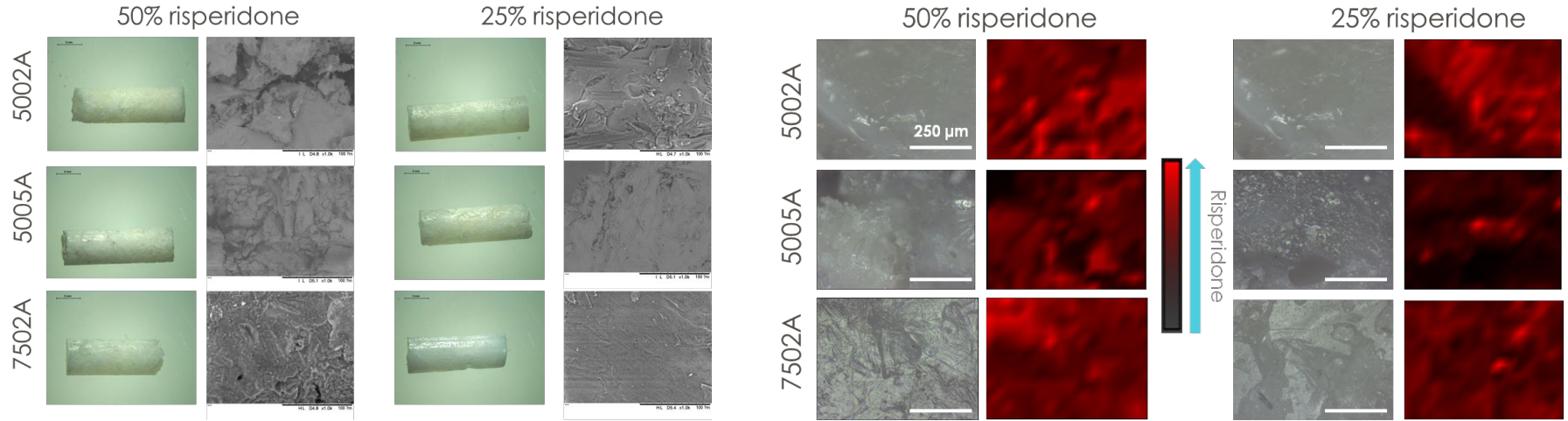




# implant characterisation – DSC + FTIR

Risperidone


↑ drug loading / ↑ drug crystallinity

L:G ratio or Mw / not a clear effect  
on drug crystallinity

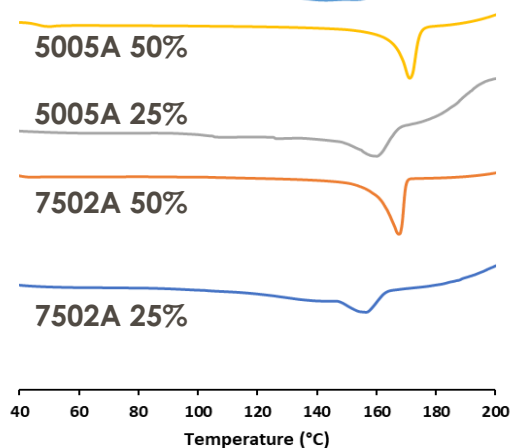

Risperidone

risperidone peaks can be clearly  
seen in the implants

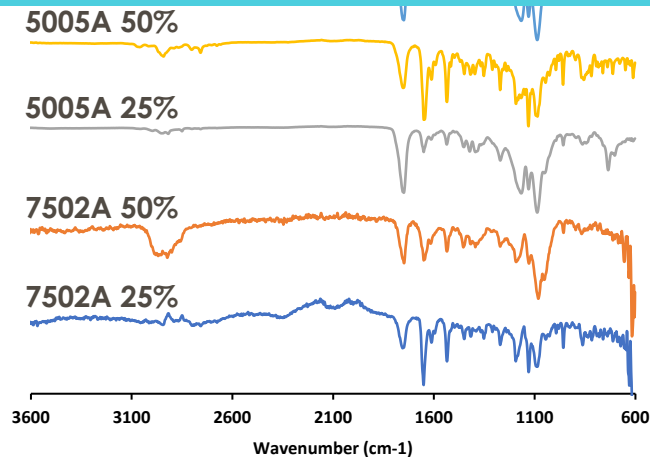





Leistritz  
EXTRUSION TECHNOLOGY



MeltPrep


Ashland™  
always solving

# implant performance testing (VCM)

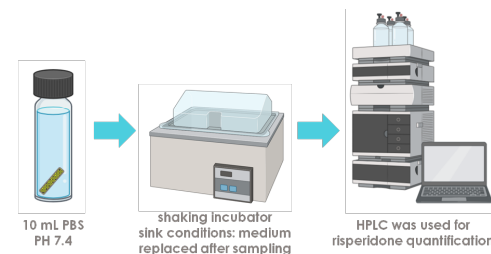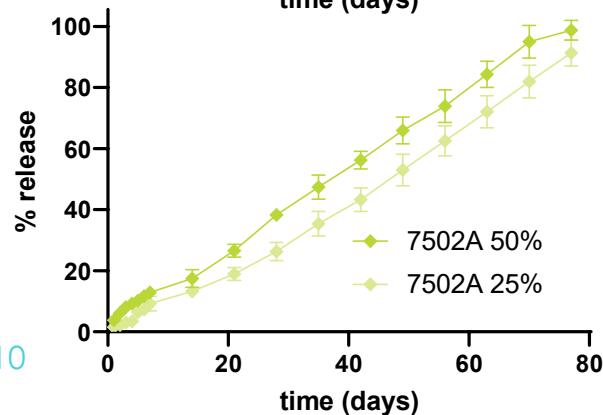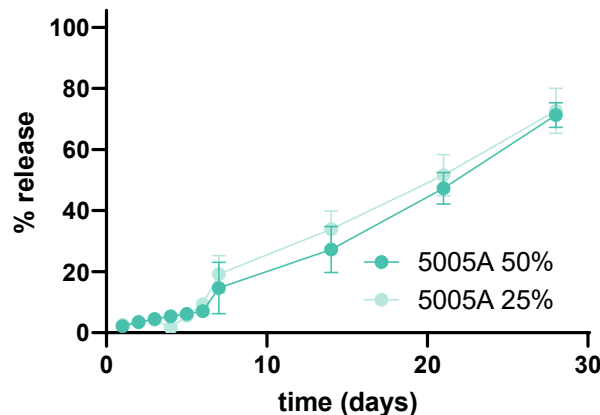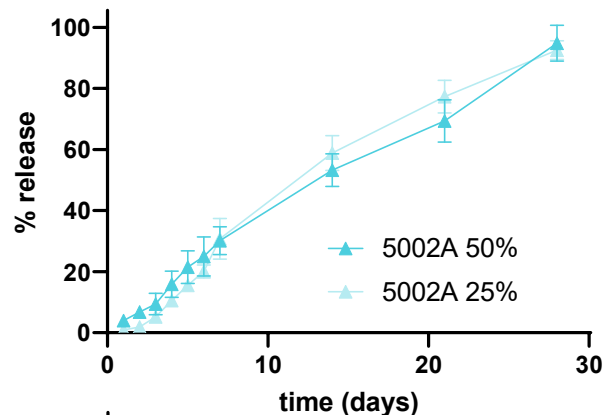

drug loading had limited effect on % release rate  
↑ drug loading / ↑ daily dose

# implant performance testing (VCM)



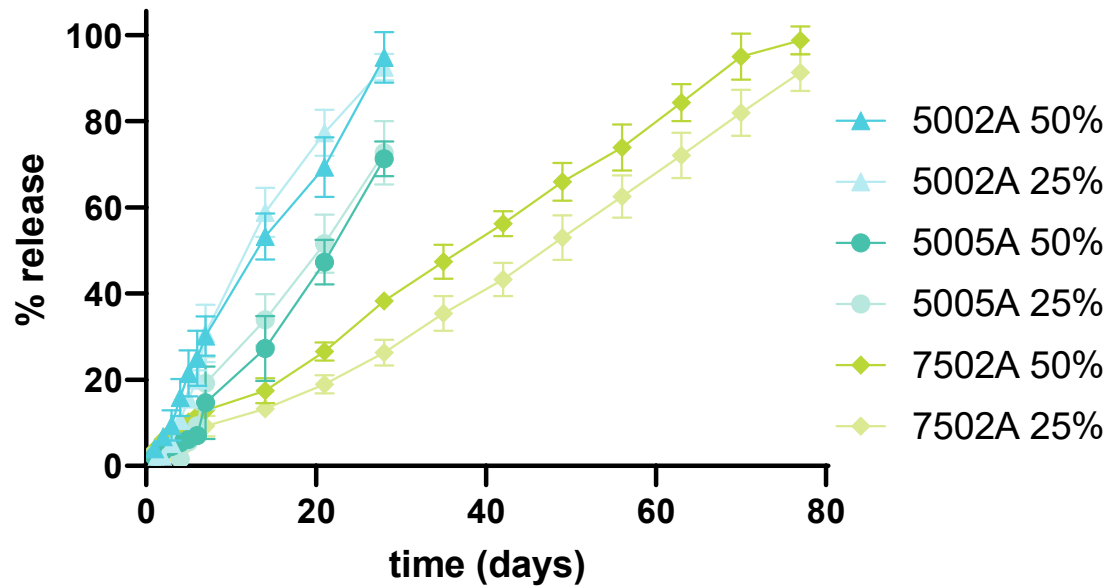


↑ Mw = ↑ release time  
(5005A slower than 5002 A)

↑ L:G ratio = ↑ release time  
(7502A slower than 5002/5A)

# scale-up via hot melt extrusion (HME)

formulations
DLG 5002 A + 50% API
DLG 5005 A + 50% API
DLG 7502 A + 50% API

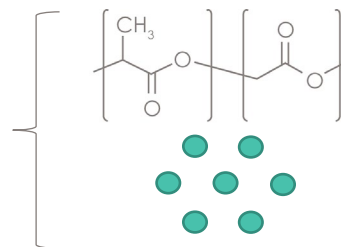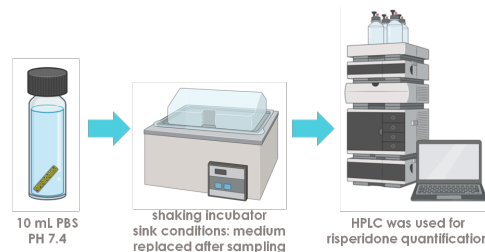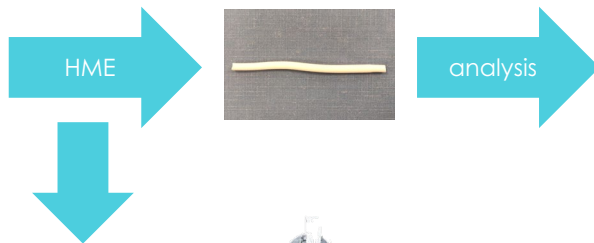

polymer

API

# implant fabrication using extruder

## Weighing

- NewClassic MS (Mettler Toledo)
- 30g polymer
- 30g API

## Mixing

- Mixomat Mini (Fuchs)
- 30min at stage 5


## Extruding

- ZSE12HPPH (Leistritz Extrusionstechnik GmbH)
- 100 g/h (gravimetric)
- 30 rpm
- 105°C at the nozzle















# implant characterisation - GPC

Sample ID	Mn (kDa)	Mw (kDa)
<b>DLG 5002 A raw</b>	<b>8</b>	<b>19.2</b>
DLG 5002 A + 50% risperidone start 1	4.7	10.7
DLG 5002 A + 50% risperidone start 2	4.9	10.9
DLG 5002 A + 50% risperidone mid 1	4.3	9.5
DLG 5002 A + 50% risperidone mid 2	4.9	10.8
DLG 5002 A + 50% risperidone end 1	5.1	11
DLG 5002 A + 50% risperidone end 2	5.2	11
<b>DLG 7502 A raw</b>	<b>10.4</b>	<b>23.1</b>
DLG 7502 A + 50% risperidone start 1	8	19.3
DLG 7502 A + 50% risperidone start 2	7.7	19.1
DLG 7502 A + 50% risperidone mid 1	7.2	18.5
DLG 7502 A + 50% risperidone mid 2	7.7	18.8
DLG 7502 A + 50% risperidone end 1	7.5	16.9
DLG 7502 A + 50% risperidone end 2	7.6	18.7
<b>DLG 5005 A raw</b>	<b>17.7</b>	<b>31.7</b>
DLG 5005 A + 50% risperidone start 1	13.1	26.0
DLG 5005 A + 50% risperidone start 2	12.4	23.5
DLG 5005 A + 50% risperidone mid 1	13.2	27.00
DLG 5005 A + 50% risperidone mid 2	13.3	26.8
DLG 5005 A + 50% risperidone end 1	13.7	27.1
DLG 5005 A + 50% risperidone end 2	13.2	26.8

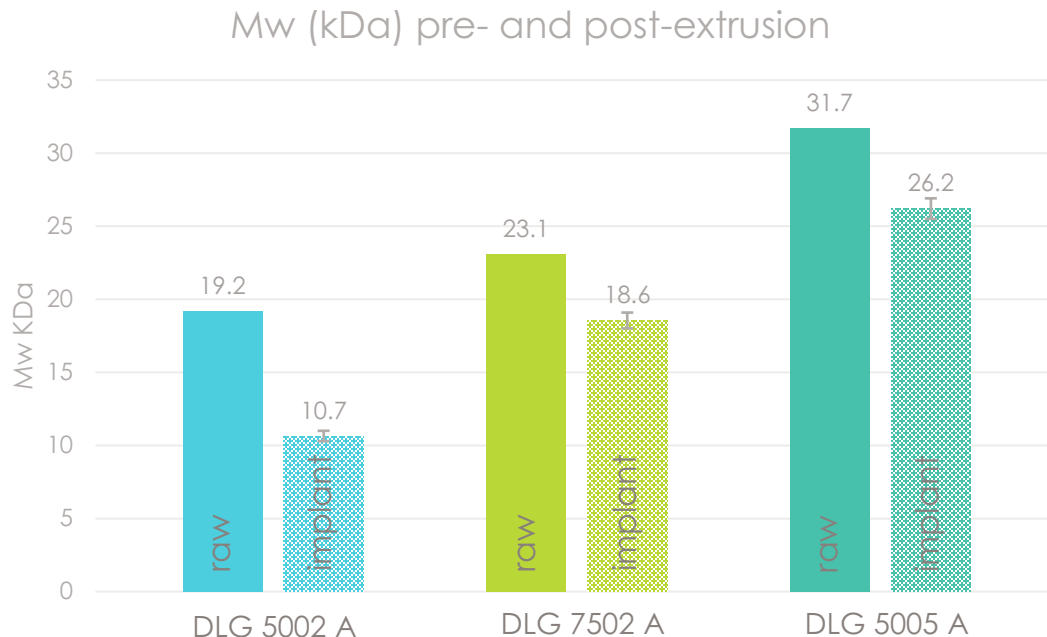

# implant characterisation – SEM & Raman


- SEM showed X
- Raman showed Y

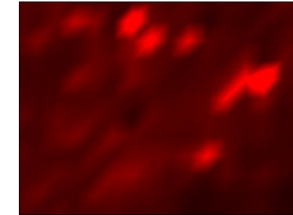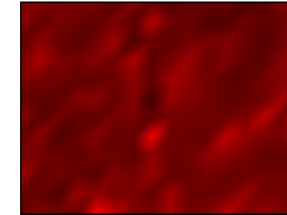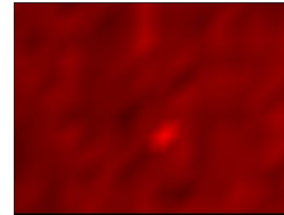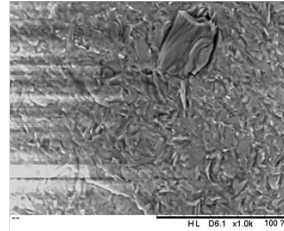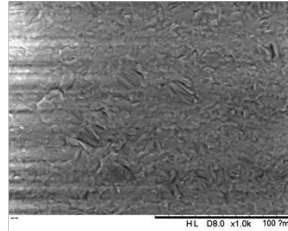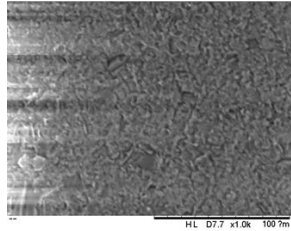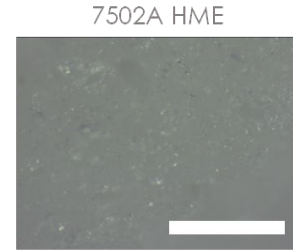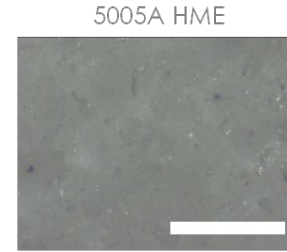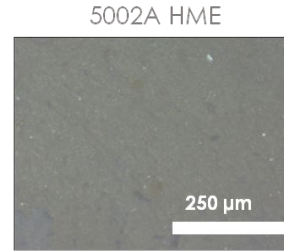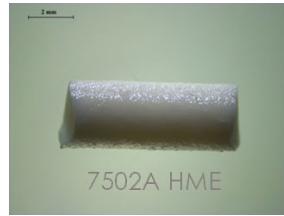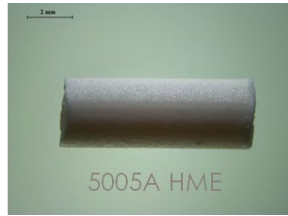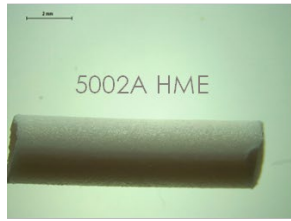


# implant comparison – FTIR & DSC

VCM and HME fabricated implants are very similar


HME

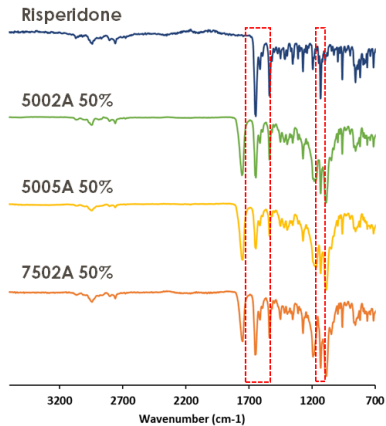


VCM

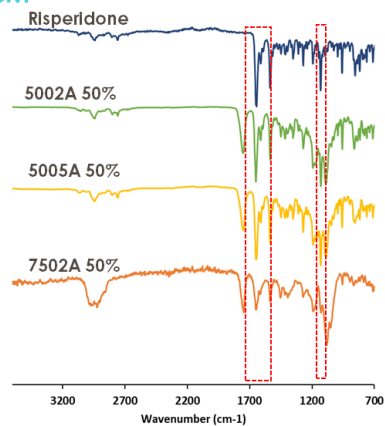

HME



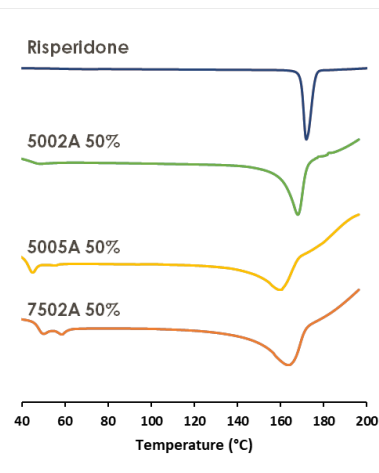


VCM

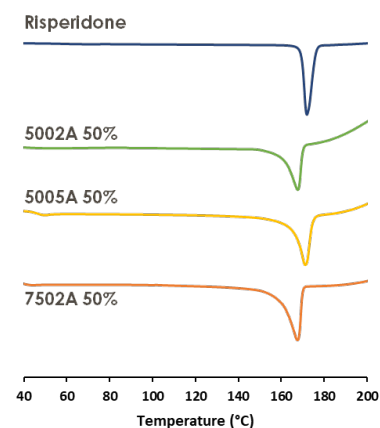



# implant performance testing

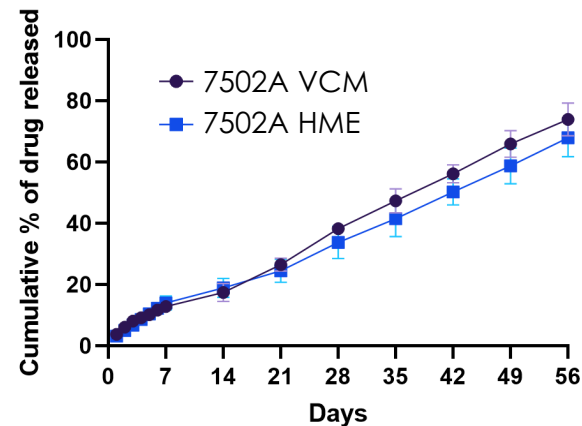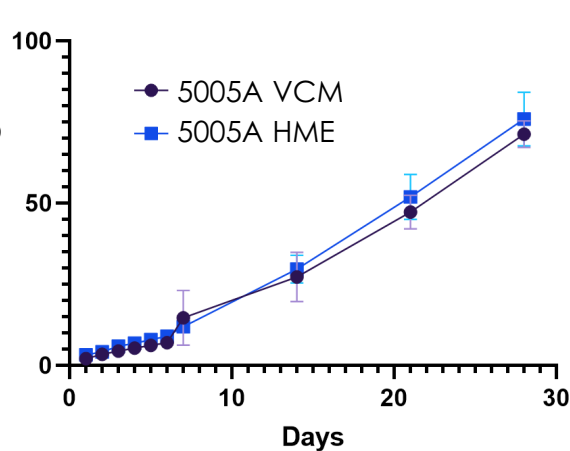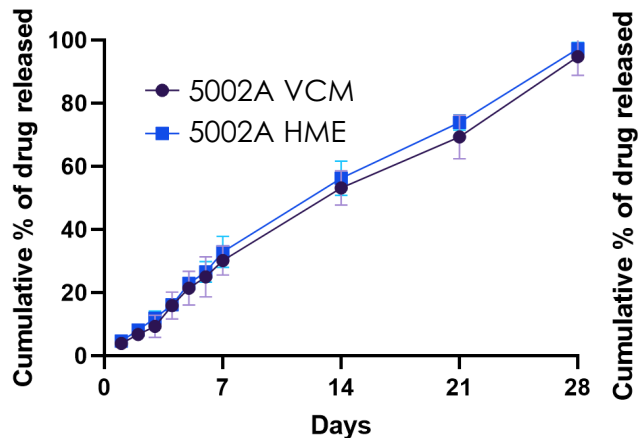

release from VCM implants = release from HME implants


# contact us



Patrick Duffy  
patrick.duffy@ashland.com



Jonas Lacina  
jlacina@leistritz.com

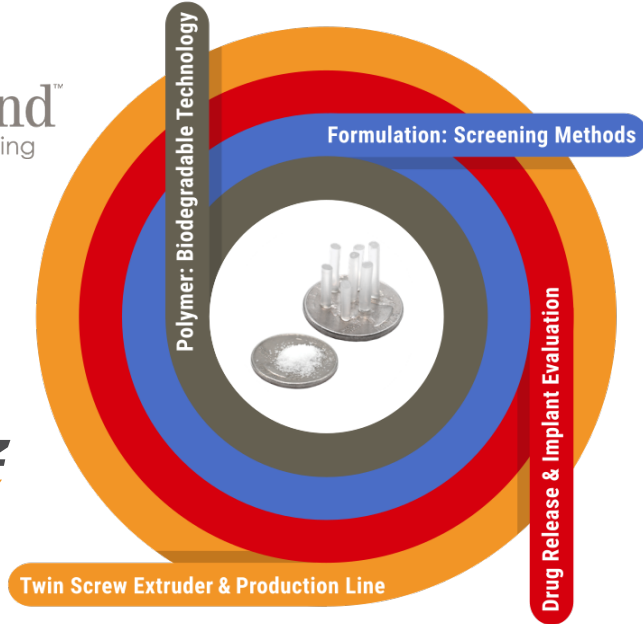

Andreas Bramböck  
andreas.bramboeck@meltprep.com



Eneko Larrañeta  
e.larraneta@qub.ac.uk



# disclaimer

---

The information contained in this presentation and the various products described are intended for use only by persons having technical skill and at their own discretion and risk after they have performed necessary technical investigations, tests and evaluations of the products and their uses. This material is for informational purposes only and describes the scientific support for the use of the products described herein as an ingredient intended to enhance performance of an end product. **Certain end uses of these products may be regulated pursuant to rules governing medical devices or other regulations governing drug uses. It is the purchaser's responsibility to determine the applicability of such regulations to its products.** While the information herein is believed to be reliable, we do not guarantee its accuracy and a purchaser must make its own determination of a product's suitability for purchaser's use, for the protection of the environment, and for the health and safety of its employees and the purchasers of its products.

Neither Ashland nor its affiliates shall be responsible for the use of this information, or of any product, method, formulation, or apparatus described in this brochure. Nothing herein waives any of Ashland's or its affiliates' conditions of sale, and no statement, information and data is to be taken as a guarantee, an express warranty, or an implied warranty of merchantability or fitness for a particular purpose, or representation, express or implied, for which Ashland and its affiliates assume legal responsibility. We also make no warranty against infringement of any patents by reason of purchaser's use of any information, product, method or apparatus described in this presentation.

The testing information (the "Testing Information") has been gratuitously provided by Ashland. The Testing Information is based on many factors beyond Ashland's control, including but not limited to, the conditions prevailing when the testing was conducted, and in some cases, is based on data generated with development samples of the Active Ingredient. Although it is intended to be accurate, ASHLAND DISCLAIMS ANY AND ALL LIABILITY, EITHER EXPRESS OR IMPLIED. The Testing Information is confidential or proprietary to Ashland, and may not, except as provided below, be disclosed to any third party. You may not make commercial use of the Testing Information or make claims with respect to your products based the Testing Information, without the written agreement with Ashland covering such use.

® Registered trademark, Ashland or its subsidiaries, registered in various countries

™ Trademark, Ashland or its subsidiaries, registered in various countries

\*third party trademark

© 2024, Ashland





Ashland<sup>TM</sup>  
always solving